

TexQ® Decontamination Protocol for Hazardous Drugs Removal

1

Remove excess liquid from the surface using a **dry wiper**.

Discard the wiper.



2

Quarter-fold a new dry wiper.



Fold in half again to make a quarter-folded wiper.

3

With nozzle on **stream**, pump the TexQ® TX650 trigger until the wiper is well dampened, but not dripping.



4

Use **unidirectional**, parallel, overlapping strokes to wipe the surface. Use one wiper for up to 40" stroke length.

Discard the wiper.



5

Wipe the surface with a sterile **70% Isopropyl Alcohol (IPA) pre-wetted wiper** using the same wiping technique described in Step 4.

Discard the wiper.



6

Repeat Steps 2 – 5 using the same wiping technique at **ninety degrees (90°)** to the previous wiping direction.